

Instructions for Mounting Flat Samples onto the Acrylic Plate

Each acrylic plate (9x9 hole pattern) is marked with a label marking Front/Top edge of the plate to define its orientation as well as the plate number. When filling the plate with samples, please follow the convention, as shown on the diagram:

- Leave two right top holes empty: they need to be used for the distance calibration and for the air blank
- Fill the rest of the plate with samples (or blanks) using Scotch tape starting from the right column and going down (see the diagram). Fold over the ends of the tape for easy sample removal.

